



**CRYSTAL TECHNOLOGY
& INDUSTRIES, INC.**

**AEOLUS™ IS-18CA/IS-18A
and IS-18C5A/IS-18B5A
Stackable CO₂/Non-CO₂
Incubator Shakers**



Company Introduction

For over 25 years, Crystal Technology & Industries has been driven to deliver more value to the global scientific community. Our approach utilizes the experience and skills of our over 40 engineers collaborating closely with top tier research institutions worldwide to design and develop products with application specific features and cutting-edge technology that provide the scientist with more innovative solutions.

The AEOLUS IS-18CA family of Stackable Incubator Shakers feature many advanced options in CO₂ Incubation, Shaking drive technology and ease of use from Crystal Technology & Industries to help scientists. All our products undergo extensive testing to ensure superior quality. That is why every piece of equipment from Crystal Industries comes with our industry leading 2-year Warranty administered by our in-house, factory trained engineers.

Our presence is global with dedicated offices/warehouses in North America, Europe and Asia. Our goal is to consistently provide total customer satisfaction felt locally in our products and services. We look forward to serving the scientific community now and the future.



BRINGING SCIENCE AND TECHNOLOGY TO LIFE!

AEOLUS™ IS-18CA/IS-18A and IS-18C5A/IS-18B5A Stackable CO₂/Non-CO₂ Incubator Shakers

The CO₂ Incubator Shaker features Temperature, Shaking, CO₂ Concentration, and Humidity controls. Its stackable design helps maximize floor space, with the capability to stack up to three units high.

INNOVATIVE DESIGN

- Three axis eccentric balance system provides stable and reliable shaking
- Terminal feedback system ensures precise shaking speed
- Brushless DC motor is highly efficient, quiet, durable, and produces negligible heat that does not affect the chamber temperature
- Combination of Proportional-Integral-Derivative control and Fuzzy control for temperature ensure temperature accuracy within $\pm 0.1^{\circ}\text{C}$
- Program mode allows user to set up to 12 segments of varying temperature and speed
- WiFi and RS485 monitoring and control allows the user to stay in touch with the instrument even outside of the lab
- Highly efficient inverter compressor refrigeration allows for temperatures as low as 20°C below room temperature with low energy consumption (minimum temperature is 4°C)
- Infrared CO₂ sensor provides accurate and reliable CO₂ detection (IS-18CA/IS-18C5A only)
- Intuitive 7" capacitive touchscreen makes functions, such as setting alarm parameters, easy to complete
- Slide rail platform makes it easy to load and unload fixtures
- Noise reduction design keeps the instrument noise level to 55 dB
- Polyurethane foam insulation provides excellent thermal insulation



CAPACITIVE TOUCH SCREEN DISPLAY

Main Interface


1. Displays the current instrument status.
2. Light button: Press to turn the light on or off.
3. Displays the current function's running status, orbit diameter (for speed), and any alarms.
4. Plus and Minus buttons: Adjusts the corresponding function's value.
5. Displays the set value and allows user to manually input the set value.
6. Run/Stop button: Runs and stops function.
7. Edit button for programmable mode: Sets up to 12 running segments of varying temperatures and speed.
8. Running mode button: Switches between "Easy" and "Program Control" mode.
9. Run/Stop all button: Runs and stops all functions.
10. Main Settings button: Enters settings page (image right), which features specific settings for the user to edit.
11. Function icon buttons: Graphically displays the real-time function data.



Control Features


- Powerful Data storage: Data is stored for more than two (2) years. Recorded in 1 minute intervals.
- Real-time data view: On the display, operator can view the running data from latest 7 days.
- Data export: Easily export more than 2 years (912 days) of parameter data via USB.
- Optional Remote WiFi / Ethernet monitoring accessory and smart phone app enabled remote monitoring and control.
- Various Alarms: Alarms can be armed to activate when parameter values deviate outside of users set bounds.
- Alarm recording: Alarm information is stored and viewable.
- Easy Calibration: Calibrating functions only require inputting the difference between set and actual values.

Overview

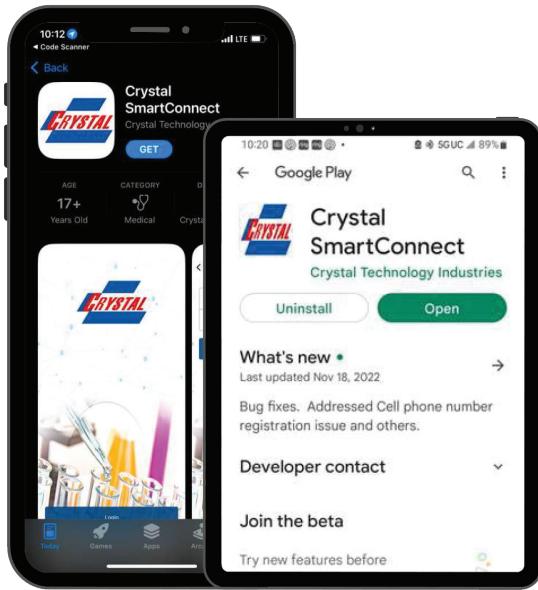

Remote Monitoring APP

Crystal Industries' Remote Monitoring accessory and app gives customers the ability to control their incubator right from their phone or tablet via WiFi Connection. Features of the incubator app include the ability to set and change the incubator's parameters (including temperature and CO2 concentration). Text notifications are sent when an alarm is activated. This unique app makes it possible to stay conveniently connected with CTI incubator and ongoing projects, whether you are in the lab or not.

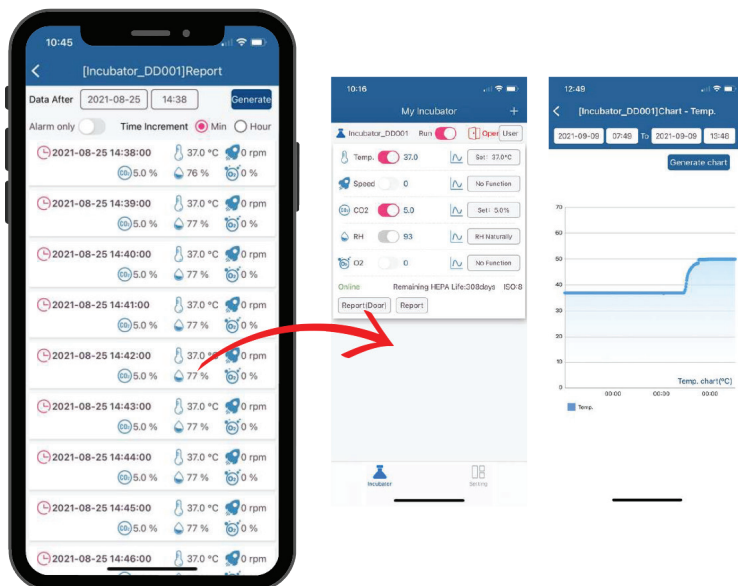

Overview

- **High Resolution 7-inch Color Touch Screen:** Modular display presents instrument status and function values in a clear and concise manner.
- **Capacitive Touch Screen:** The capacitive screen is more responsive and durable than a resistance screen as there is no reaction delay when the screen is pressed.
- **Pull Out Rail Platform:** Operators can easily pull the platform out to load or unload samples. Only one locking handle in the middle of the platform makes the process fast and convenient. Universal Platform is shipped pre-installed in the unit.
- **Excellent Temperature Control Performance:** When the incubator is at a stable temperature of 37° C, the uniformity is $\pm 0.1^{\circ}$ C.
- **Strong Cooling Capacity:** The incubator chamber can be cooled to 15° C below the ambient temperature (Lowest temp is 4° C).
- **Inverter Compressor:** Makes cooling a quiet, durable, and highly efficient process while also saving energy.
- **Three Units Can Be Stacked:** This product can be purchased as a single, double, or triple stack.
- **Bi-Directional Ultraviolet Sterilization:** The chamber is equipped with a set of UV sterilization light components to enhance the germicidal irradiation effect and reach all areas of the chamber.
- **Ultra Quiet Design:** The noise level is less than 55dB when incubator is running and cooling (loudest process), providing a quiet environment for the operator.
- **Slow Acceleration and Deceleration:** Shaking function starts and stops slowly to prevent large shear forces on the bacteria or cells.
- **Efficient, Durable, and Lightweight Insulation:** Polyurethane foam insulation provides advanced thermal insulation. The material is also air-tight, waterproof, lightweight, shockproof, electrically insulated, heat resistant, cold resistant, and solvent resistant.

AEOLUS™ CO2/Non-CO2 Incubator Shakers


AEOLUS™ IS-18CA/IS-18A CO2/Non-CO2 Incubator Shaker

The CO2 Incubator Shaker features Temperature, Shaking, CO2 Concentration, and Humidity controls. Its stackable design helps maximize floor space, with the capability to stack up to three units high.





◆ Features

- 7" High-Res Color Touch Screen with intuitive graphic interface
- Chamber design features smooth rounded internal corners for easy cleaning
- Capacity is 6 x 5L Thompson shaker flasks
- Infra-Red(IR) CO2 sensor for precise and reliable CO2 detection
- Optional WiFi and RS485 monitoring and control
- Audible and visual alarms
- Mirror finish, pharmaceutical-grade stainless steel chamber
- Stackable design up to 3 units
- Choose your Custom Stack, choice of CO2 and NonCO2, 25 or 50mm Orbit
- 25mm and 50mm Orbit Models available
- 2 Year Warranty
- Units shipped pre-stacked on single pallet, single, double or triple stack (Stand sold separately)

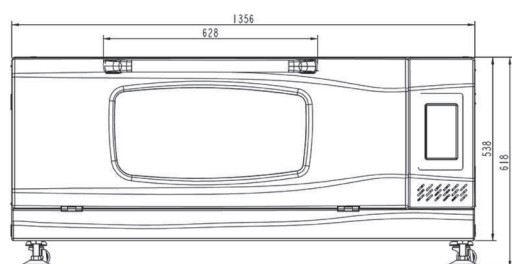


Single Unit Dimensions
Front View

Single Unit Dimensions
Side View

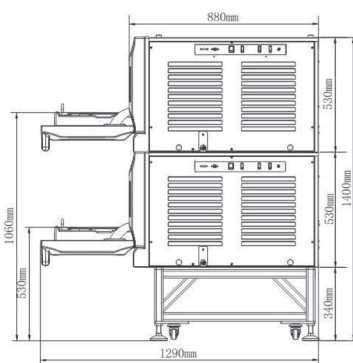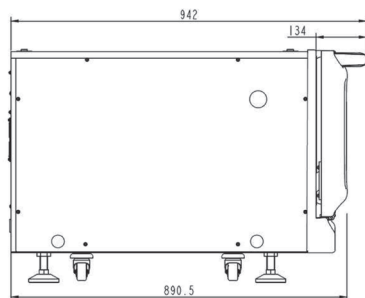

Double Stack with IS-A118 Stand
Dimensions

Triple Stack with Down-Swing Doors
Dimensions

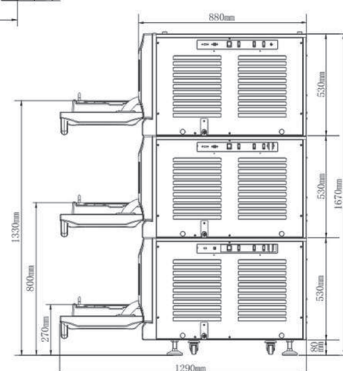

AEOLUS™ CO2/Non-CO2 Incubator Shakers

PRODUCT SPECIFICATIONS

Model	IS-18A	IS-18B5A	IS-18CA	IS-18C5A
Product Name	AEOLUS™ Incubator Shaker	AEOLUS™ Incubator Shaker (50mm orbit)	AEOLUS™ CO2 Incubator Shaker	AEOLUS™ CO2 Incubator Shaker (50mm orbit)
Platform Dimensions (WxD)	900 x 560 mm (35.4 x 22 in)			
Max Capacity*	Clamp capacity: 25mLx134 / 50mLx134 / 100mLx60 / 150mLx60 / 200mLx60 / 250mLx60 / 500mLx32 / 1000mLx22 / 2000mLx15 / 5000mL (Shaker Flask) x 6 Capacity with Sticky Mats: 25mLx187 / 50mLx187 / 100mLx112 / 150mLx104 / 200mLx84 / 250mLx76 / 500mLx40 / 1000mLx28 / 2000mLx15 / 5000mL (Shaker Flask) x 6			
Capacity	302 L			
Max Load	20 kg (44.1 lb)			
Temperature Range	Room Temp. Minus 20°C~60°C (Lowest Temp is 4°C)			
Temperature Accuracy	±0.1°C (@37°C)			
Temperature Uniformity	±0.1°C (@37°C)			
Temperature Stability	±0.1°C (@37°C)			
Temp Recover Time After Door is Opened for 30s	≤6min (@37°C)			
Speed**	0 (stop), 30~300 RPM			
Speed Accuracy	±1 RPM			
Orbit Diameter	Ø25mm (Ø1.0in)	Ø50mm (Ø2 in)	Ø25mm (Ø1.0in)	Ø50mm (Ø2 in)
CO2 Sensor	NA		Infrared	
CO2 Control Range	NA		0~20%	
CO2 Control Accuracy	NA		+0.1%	
CO2 Recover Time After Door is Opened for 30s	NA		≤6min (@5%)	
Relative Humidity Type	NA		Active	
Max Humidity	NA		95%	
Programmable Segments	12 Segments			
Timer	0-999hr 59min			
Illumination	LED Lamps on Left and Right Walls			
Disinfection	UV Lamps on Left and Right Walls			
Door Opening Direction	Down-Swing (Optional up-swing door for top unit of triple stack)			
Internal Working Dimensions (WxDxH)	900 x 560 x 308mm (35.43 x 22.05 x 12.13in)			
External Dimensions (WxDxH)	1355 x 855 x 535 (+80mm w/ feet) mm (53.35 x 33.66 x 21.06 (+ 3.1in w/ feet) in)			
Internal Walking Height	308mm (12.1in)			
Net Weight	195 kg (430 lb)			
Display	7" Color Capacitive Touch Screen			
Parameter Protection	Password Protected			
Data Stored	Time, Temperature, Speed, Instrument Status		Time, Temperature, Speed, CO2 Concentration, Humidity, Instrument Status	
Data Storage Interval	1 Minute			
Screen Data Display	Displays Most Recent 7 Days of Data			
USB Data Export	Most Recent 912 Days of Data			
Automatic Defrost	Yes			
Noise Level	≤55dB (When Compressor is Running)			
Auto-Recovery after Power Interruption	Yes			
Power supply	AC 110V±11V, 60Hz			
Power when Heating	≤800w			
Power when Cooling	≤300w			
Max Power	≤1000w			
Mobile Monitoring and Control by WiFi	Optional through IS-A64W Accessory and App			
Warranty	2 Years Parts and Service			
Certification	CE			

*Sticky Mat can be cut to fit platform

**Note: Double Stack units is recommended that the Shaker Speed of the top unit <250 rpm

Triple Stack units is recommended that the Shaker Speed of the middle and top unit < 250 rpm

AEOLUS™ CO2/Non-CO2 Incubator Shakers

Accessories



Test Tube Rack



Adjustable Test Tube Rack



Mini Universal Spring Platform



Infusion Bottle Clamp



Bioreactor Tube Holder



Flask Clamp



Stand for IS-18CA/IS-18A Incubator Shakers (Single or Double Stack)



96-Well Microplate Holder



Sticky Mat



CO2 Cylinder Pressure Regulator



Water Carboy



Multi-Microplate Holder



Fixture for Separating Funnel



Visor Cover

Item Number	Item Type	Specifications	Max Capacity
IS-A50	Flask Clamp	Holds 25mL Flask	134
IS-A1		Holds 50mL Flask	134
IS-A2		Holds 100mL Flask	60
IS-A3		Holds 150mL Flask	60
IS-A4		Holds 200mL Flask	60
IS-A5		Holds 250mL Flask	60
IS-A6		Holds 500mL Flask	32
IS-A7		Holds 1000mL Flask	22
IS-A9		Holds 3000mL Flask	12
IS-A10		Holds 5000mL Flask	6
IS-A51	Tube Rack	Holds 50 Tubes x 1.5mL/2.0mL/5mL	24
IS-A11		Holds 40 Tubes x Ø14mm	9
IS-A12		Holds 40 Tubes x Ø16mm	9
IS-A23		Holds 40 Tubes x Ø18mm	9
IS-A24		Holds 27 Tubes x Ø22mm	9
IS-A42		Holds 21 Tubes x Ø30mm	9
IS-A20	Adjustable Tube Rack	Holds 40 Tubes x Ø14mm	9
IS-A21		Holds 40 Tubes x Ø16mm	9
IS-A30		Holds 40 Tubes x Ø18mm	9
IS-A31		Holds 27 Tubes x Ø22mm	9
IS-A32		Holds 21 Tubes x Ø30mm	9
IS-A37		Holds 24 Tubes x Ø30mm	6
IS-A33	Infusion Bottle Clamp	Holds 500 ml Infusion Bottle	40
IS-A34		Holds 1000 ml Infusion Bottle	28
IS-A13	Microplate Holder	96-Well Microplate Holder	30
IS-A175	Multi-Microplate Holder	Holds up to 6 Standard Microplates	30
IS-A167	Deep Well Plate Holder	Holds 2 Deep Well Plates	30

Item Number	Item Type	Specifications	Max Capacity
IS-A35	Fixture holds Separating Funnel	Holds Separating Funnels of ±500ml	4
IS-A47	Sticky Mat	140 x 140mm (5.51 x 5.51in)	24
IS-A122		280 x 280mm (11.02 x 11.02in)	6
IS-A63	Spring Frame	900 x 540 x 80mm	1
IS-A27		350 x 240 x 80mm	4
IS-A22		450 x 400 x 80mm	1
IS-A14		428 x 295 x 80mm	2
IS-A90	Water Carboy	For IS-18CA humidity function; Holds 15L; 1 can serve up to 3 incubator shakers	1
IS-A64W	Remote Monitoring and Control	Connects each unit to users' smart phones via a WiFi or wired internet connection and an app available on the Apple Store and Google Play	1
CIA25	CO2 Cylinder Pressure Regulator	Max Inlet Pressure: 3000 psi Max Outlet Pressure: 60 psi	—
IS-A118	Stand	Adds 34mm (1.339in) of height Recommended for use with Single and Double Stacks	—
IS-A166	Bioreactor Tube Holder	Holds 450ml Bioreactor Conical Tubes	6
IS-A178		Holds 600ml Bioreactor Conical Tubes	6
IS-A176	Visor Cover	Blocks ambient light from entering chamber	1



Crystal Industries Sales & Product Support:

1600 S. Airport Dr, Suite 100, McKinney, TX 75069

972-934-2525

crystalindustries.com

